





AQ400

DISCRETE ANALYZER FOR ENVIRONMENTAL TESTING

AQ400 is a flexible analyzer that uses the principle of discrete analysis where each test occurs in a separate or discrete reaction vessel.



AQ400 is ideal when many and varied tests are needed on different samples and/or individual results are needed immediately.

USEPA, ASTM, ISO

and other international regulatory compliant methods are available.

METHODS INCLUDE

Alkalinity

Ammonia

Chloride

Cyanides

Nitrate/Nitrite

Nitrite

Phenol

Phosphate, ortho


Phosphorus, total

Silicate

Sulfate

Total Kjeldahl Nitrogen

PLUS MANY MORE

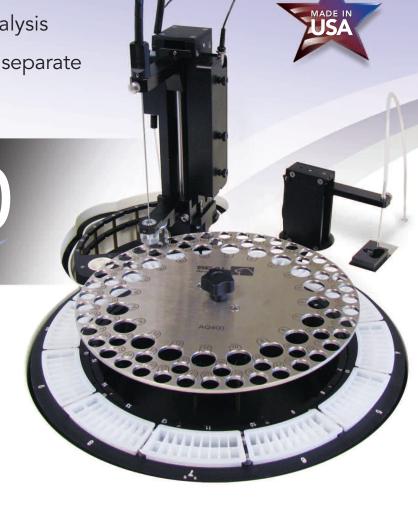

HOW DOES THE AQ400 WORK?

The AQ400 robotic sampling arm works in conjunction with a stepper motor-driven syringe that is responsible for aspirating, dispensing and mixing accurate and precise quantities of sample and reagent in miniaturized test tubes called reaction wells.

The sample and reagents are incubated in heated reaction wells until reaction is complete. A single aliquot is then transferred into a 10mm pathlength optical glass cuvette. The absorbance is read on the stationary reactant to ensure the best possible signal to noise ratio.

Each sample is read is the same position in front of the detector in the 10mm glass cuvette, similar to SEAL colorimetric flow systems, known for their high reproducibility and lowest detection limits. This eliminates issues of reaction well variability and scratching found in direct read discrete systems so improves reproducibility. In the SEAL AQ series, liquid is moved not the tray - fewer moving parts equals higher reliability.

Once the absorbance is read, the glass cuvette is thoroughly cleaned and checked, ensuring no carryover or cross contamination.

As a market leader, SEAL has over 1,000 applications available and under continual development. Markets include water, wastewater, soil, plant, fertilizer, food and beverage. Please contact us for your specific application.

AQ400 – The latest in discrete analyzer technology Designed specifically for the environmental market

ADVANTAGES AND BENEFITS



- ~ 100% optical glass stop-flow cuvette
- ~ 10 mm optimum path length
- ~ Different path lengths available



- Low cost, disposable wells used for each discrete reaction
- Constant heating and programmable reaction time ensure reaction reaches completion
- Automatic reagent level sensing verifies sufficient reagent volume



- Optional integrated cadmium coil reduction for nitrate/nitrite determination
- Cadmium coil is sealed and valve controlled
- In situ coil regeneration is fully automated



- Unique probe washer for cleaning sample probe between sample and reagent
- ~ Eliminates cross contamination
- ~ Ideal for wastewater



- Highly flexible software designed with user input
- ~ QCPro™ Data Quality System – allows the user to specify QC types, limits and corrective actions

- True unattended operation including ability to run overnight
- Automated standard preparation and dilution of over range samples
- Tests programmable per sample to reduce analysis time
- Add samples after a run has started
- Total volume per test only 500 600 μL
- Different size sample trays are available to accommodate different workloads
- Segregated chemical waste and wash minimizes waste disposal
- LIMS compatible export in .csv format



The AQ400 can be used as a standalone spectrophotometer.

The optional vial adaptor can accept vials of different sizes for reading tests such as COD.



www.seal-analytical.com

Colorimetric Nutrient Analyzers

DISCRETE ANALYZERS







AQ270

AQ300

AQ400

SEGMENTED FLOW ANALYZERS







QuAAtro39

AA3

AA500

50 Years' Experience in Environmental **Analysis Built into Every Analyzer**

50 years' experience in designing, developing and manufacturing automated wet chemistry analyzers specifically for very low detection levels in environmental applications has helped SEAL to apply the most useful, easy to use features into the SEAL range of Discrete and Segmented Flow analyzers. The SEAL analyzers are widely acknowledged as the best for environmental analysis, giving you everything you need to achieve equal or superior results to the manual and approved laboratory methods the SEAL analyzer replaces.

Digestion Systems



BD50

FOR METALS AND TKN. **TP DIGESTION**



SmartBlock II



DEENA 3

SEAL Analyzers are monitoring environmental samples in every corner of the globe. They are manufacturered in the USA, Germany and the Netherlands. Engineering and chemistry support is provided from SEAL global facilities in USA, Germany, England, the Netherlands and China along with a worldwide network of specialist distributors.

COMPREHENSIVE SUPPORT

We offer comprehensive applications, technical service and software support.

INCLUDING

- A choice of preventative maintenance and service contracts to meet your specific requirements
- In-house and online training
- Guaranteed availability of genuine consumables and spare parts
- Adaptation of methods to specific requirements such as matrix, range or detection limit
- Continuous in-house development of software to incorporate new customer requested features

Robotic Handling Systems

SEAL Robotic Minilab systems for automating sample (pre) treatment in the laboratory improving your sample handling efficiency. Typical applications include BOD, pH, COD, Alkalinity, and conductivity measurements with options such as decapping/capping, sample splitting, and filtration. Call us about your laboratory needs and we will design a robot to suit you.





www.seal-analytical.com

SEAL Analytical is a global company with offices worldwide - contact us at:

SEAL Analytical, Inc.

sales@seal-us.com

6501 West Donges Bay Road Mequon, WI 53092 **United States** Tel: +1 (262) 241 7900 Fax: +1 (262) 241 7970

SEAL Analytical Ltd.

3 Talisman Business Centre Duncan Road Park Gate Southampton, S031 7GA United Kingdom Tel: +44 (0) 1489 864 400 sales.uk@seal-analytical.com

SEAL Analytical Netherlands

ROHASYS BV Provincienbaan 4 5121 DL Rijen The Netherlands Tel: +31 161 240152 Fax: +31 161 240153 info@rohasys.com

SEAL Analytical GmbH

Werkstrasse 5 D-22844 Norderstedt Germany

Tel: +49 (0)40 60 9292 9-00 Fax: +49 (0)40 60 9292 9-02 info.germany@seal-analytical.com

SEAL Analytical Shanghai

Room 413, 12th Building, No. 128 Xiangyin Road, Shanghai, 200433 China

Tel: +86 21 3362 5002 Fax: +86 21 3362 5002